Technical Working Party on Testing Methods and Techniques

TWM/2/17

Original: English Date: April 8, 2024

GUIDELINES FOR THE VALIDATION OF A NEW CHARACTERISTIC-SPECIFIC MOLECULAR MARKER PROTOCOL AS AN ALTERNATIVE METHOD FOR OBSERVATION

Document prepared by experts from France, Italy and the Kingdom of the Netherlands

Disclaimer: this document does not represent UPOV policies or guidance

Table of Contents:

ASS	SOCIATED DOCUMENTS	2
I.	OBJECTIVES OF THESE GUIDELINES	2
II.	SCOPE OF THESE GUIDELINES	2
III.	PERFORMANCE CRITERIA FOR A NEW MOLECULAR MARKER BASED PROTOCOL	2
S	Specificity	2
	Definition Requirement	2 2
S	Sensitivity and limit of detection	3
	Definition Requirement How to evaluate it?	
F	Repeatability	3
	Definition (based on ISO 16 577:2016) Requirement How to evaluate it?	
F	Reproducibility	3
	Definition (based on ISO 16 577:2016) Requirement How to evaluate it?	
F	Robustness	4
	Definition Requirement How to evaluate it?	
IV.	VALIDATION REPORT	4
V. PRC	STANDARD PROTOCOL FOR CHARACTERISTIC-SPECIFIC MOLECULAR MARKER DTOCOL	5
VI.	FOLLOW-UP SURVEY AFTER APPROVAL	8
ANN	NEX Presentation by CREA, GEVES and Naktuinbouw	

ASSOCIATED DOCUMENTS

- 1. Please note that the parts highlighted indicate text quoted from the documents below:
 - TG/1/3: General Introduction to the Examination of Distinctness, Uniformity and Stability and the Development of Harmonized Descriptions of new Varieties of Plants
 - TG/44: Guidelines for the conduct of tests for distinctness, uniformity and stability for Tomato
 TGP/12: Guidance on Certain Physiological Characteristics
 - TGP/15: Guidance on the Use of Biochemical and Molecular Markers in the Examination of Distinctness, Uniformity and Stability (DUS)
 - UPOV/INF/17 Guidelines for DNA-Profiling: Molecular Marker Selection and Database Construction
 - UPOV/INF/18 Possible use of Molecular Markers in the Examination of Distinctness, Uniformity and Stability (DUS)
 - TWV/54/7 + Add Use of molecular techniques in DUS examination

I. OBJECTIVES OF THESE GUIDELINES

2. The purpose of these guidelines is to elaborate the principles contained in the General Introduction (document TG/1/3), and its associated TGP documents, into detailed practical guidance for the harmonized validation of a new method based on molecular marker before its use as an alternative test. Performance criteria required for the validation are described and guidance on their assessment is given. These guidelines also describe a standard protocol with mandatory and optional chapters. Survey after acceptance is also described.

3. If a different technique is used, the laboratory must validate its method in comparison to the reference method (to show that the alternative technique gives the same results).

II. SCOPE OF THESE GUIDELINES

All crops

Characteristic-Specific Molecular Markers For the examination of Distinctness, Uniformity, and Stability (DUS).

III. PERFORMANCE CRITERIA FOR A NEW MOLECULAR MARKER BASED PROTOCOL

Specificity

Definition

4. Correlation between the genotype and the phenotype, *i.e.* reliability of the link between the marker and the characteristic.

Requirement

5. In principle 100% of correlation between the genotype and the phenotype. If the correlation is less than 100% a follow-up test(s) should be performed to ensure the reliability of the results. A decision rule can be used in that case. Less than 100% correlation can be caused by other genetics. It can also suggest that the non-correlation is caused not by the marker but by external factors in the phenotypical observations (*e.g.* biotest for a disease resistance).

How to evaluate it

6. Number of varieties: "To start the marker selection process an appropriate number of varieties (development set) is needed to reflect at the most the diversity observed within the group/crop/species/type for which the markers are intended to be discriminative."

7. Varieties should represent the different states of expression (if known varieties with heterozygous and homozygous state), coming from different seed companies, with different genetic background of the characteristic and different types. Well phenotypically characterized varieties for the trait of interest should be used when available.

8. Number of plants per variety: At least one plant per variety if there is available well phenotypically characterized varieties. If not, the number of plants must be the same as for the morphological observation described in the UPOV guideline.

9. The specificity can be assessed within one laboratory.

Sensitivity and limit of detection

Definition

10. The limit of detection is defined as the minimal quantity of the target that can be reliably detected.

11. In case of analyses performed on bulk samples (*e.g.* pool of different plants of the same variety) the sensitivity is critical and must be assessed. For the use on individual plants, the quantity of the target is not critical and this performance criterium is optional.

Requirement

12. In the case of the pool, the requirement would be to detect at least one off-type in the pool.

How to evaluate it?

13. To use artificial samples by mixing one off-type to a pool to check the sensitivity of the detection.

Repeatability

Definition (based on ISO 16 577:2016)

14. *"Repeatability; where identical test results are obtained with the same method, on identical test items, in the same laboratory, by the same operator, using the same equipment within short intervals of time."*

15. For qualitative methods, accordance is equivalent to the repeatability of quantitative methods (Langton *et al.*, 2002).

Requirement

16. Ideally 100%, a performance \geq 90% is generally accepted. If the repeatability of the reference method is published the repeatability of the alternative method should be at least equivalent.

How to evaluate it?

17. The repeatability can be evaluated within one laboratory.

18. At least three technical replicates drawn from a same plant (three independent DNA extractions). To include at least all expected types of genotype.

Reproducibility

Definition (based on ISO 16 577:2016)

19. *"Reproducibility; where test results are obtained with the same method, on identical test items, within the same laboratory or between different laboratories, with different operators, using different equipment*" at different times.

20. For qualitative methods, concordance is equivalent to the reproducibility of quantitative methods (Langton *et al.*, 2002).

Requirement

21. Ideally 100%, a performance \geq 90% is generally accepted. If the reproducibility of the reference method is published the reproducibility of the alternative method should be at least equivalent.

How to evaluate it?

22. Reproducibility should be assessed between different laboratories by an interlaboratory validation study (Ring-test) with coded samples of known genotypes. All expected types of genotype should be included.

23. The ring-test should involve at least, three different laboratories including at least two different examination offices (*e.g.* in the INVITE project 3 examination offices were involved in the validation test). If possible, experienced laboratories familiar with the species and the technique should be involved. If not, a training can be organized ahead of the ring-test with un-coded samples. Laboratories can participate in a ring-test on voluntary basis. In case there are no volunteers, the reproducibility can be determined within one laboratory.

24. All laboratories must follow the protocol to be validated. In the protocol compulsory and optional parts can be defined by the validation team. For example, see the protocol CPVO/TP-044/4-Rev. where compulsory and optional steps were defined.

25. Number of varieties: To include at least all expected types of genotype.

26. Guidelines/Norms on interlaboratory studies can be followed: ISO 13495 Foodstuffs - Principles of selection and criteria of validation for varietal identification methods using specific nucleic acid, ISO 17043 Conformity assessment - General requirements for proficiency testing, EPPO pm7-122-2 Guidelines for the organization of interlaboratory comparisons by plant pest diagnostic laboratories, ISTA TCOM-P-10-Validation of seed health methods and organization and analysis of interlaboratory comparative tests (CT)... The validation team can cite the followed guidelines in its report.

<u>Robustness</u>

Definition

27. "Robustness; a measure of its capacity to remain unaffected by small, but deliberate deviations from the experimental conditions described in the procedure parameters and provides an indication of its reliability during normal usage" (e.g. change of DNA extraction method or change of real time machine).

Requirement

28. Ideally 100%, if less that means that the method is not robust to a change of one parameter and this should be indicated in the protocol as a mandatory step (*e.g.* a change of a mastermix that would be critical).

How to evaluate it?

29. It is optional to assess, and robustness is evaluated partially during the ring test (reproducibility), (different laboratories, equipment, machinery, persons, etc.).

IV. VALIDATION REPORT

30. The validation report and results must be peer-reviewed by two (preferably 3 if the reproducibility was done within one laboratory) of the examination offices. Reviewing is on voluntary basis but preferably perform review by laboratory familiar with the species and the method.

31. During the reviewing process, the reviewers can require extra validation data in concertation with the validation team.

Content of the validation report

- Raw data generated during the different steps of the validation process
- Detail protocol with optional and compulsory steps defined
- Performance criteria assessment
- Conclusion

Publicity

32. The validation report should be available upon request. In the new protocol the validation process should be mentioned with the contact examination office. In some particular cases, *e.g.* a "trade secret protocol" (cytoplasmic male sterility in cabbage), the protocol and the validation report could not be shared outside of the examination offices.

V. STANDARD PROTOCOL FOR CHARACTERISTIC-SPECIFIC MOLECULAR MARKER PROTOCOL

33. Compulsory elements are indicated in the column "essential information", the other elements may be used depending on the characteristic test protocol. If a laboratory wants to adapt/modify/change a mandatory chapter or element of a mandatory chapter it must validate its method in comparison to the reference method (to show that you obtain the same results as the published method).

Table 1: Standard characteristic-specific molecular	marker protocol	(see document	TWV/54/7 '	Use of
molecular techniques in DUS examination	n". Modifications	are highlighted	in grey)	

Chapter	Elements in a Standard characteristic- specific molecular marker protocol	Example	Essential information for harmoni- zation	Remark
1	characteristic	Resistance to Tomato mosaic virus (ToMV) See TG/44/11/rev3 – Ad 51: ii DNA marker test	YES	
2	Genes and alleles	See TG/44/11/rev3 – Ad 51: ii DNA marker test add 2	YES	Need to avoid dominant marker or presence/absence marker otherwise the robustness should be assessed
	Targeted gene(s)	Resistance Gene Tm2		a) file(s) containing the DNA sequence information (order of nucleotides)
		Arens, P. et al (2010)		b) reference to DNA information in public databases (like GeneBank)
2.1			YES	c) reference to (scientific) publications in which the DNA sequence information of the states of expression of the characteristic is revealed.
				d) reference to a particular position on the published reference genome version.
		Tm2 and Tm2 ²		a) file(s) containing the DNA sequence information (order of nucleotides)
		Arens, P. et al (2010)		b) reference to DNA information in public databases (like GeneBank)
2.2	Allele corresponding to state 1		YES	c) reference to (scientific) publications in which the DNA sequence information of the states of expression of the characteristic is revealed.
				d) reference to a particular position on the published reference genome version in combination with the SNP or INDEL that is responsible for the state of expression.

Chapter	Elements in a Standard characteristic- specific molecular marker protocol	Example	Essential information for harmoni- zation	Remark
2.3	Allele corresponding to expression state n	tm2 Arens, P. et al (2010)	YES	 a) file(s) containing the DNA sequence information (order of nucleotides) b) reference to DNA information in public databases (like GeneBank) c) reference to (scientific) publications in which the DNA sequence information of the states of expression of the characteristic is revealed. d) reference to a particular position on the published reference genome version in combination with the SNP or INDEL that is
3	Primers (and probes)	See TG/44/11/rev3 – Ad 51: ii DNA marker test add 3, 3.1 and 3.2	YES	responsible for the state of expression. Primer and probe sequences, reference to accessions and sequences in public databases (Genebank numbers), literature
3.1	Primers (and probes) to detect allele '9'		YES	Primer Sequences corresponding to allele(s) for expression '9' (resistance)
3.2	Primers (and probes) to detect allele '1'		YES	Primer Sequences corresponding to allele(s) for expression '1' (susceptibility)
3.3	Primers (and probes) to detect allele 'x'		YES	Primer Sequences corresponding to allele(s) for expression 'x'
4	Format of the test			
4.1	Number of plants per genotype	>=20	YES	A minimal number of individual plants required (see 5.2.1a) the test for the marker is conducted on the same number of individual plants, with the same criteria for distinctness, uniformity and stability as for the examination of the characteristic by an observation assay (TGP 15)
4.2	Control varieties	See TG/44/11/rev3 – Ad 51: ii DNA marker test add 4.2	YES	Control varieties (same as in observation assay) as standards representing all relevant combination of alleles. For example homozygous for Allele corresponding to expression state 9 (present), homozygous for allele corresponding to expression state 1 (susceptible) and heterozygous (both alleles are present in a diploid) corresponding to either resistant, susceptible or intermediate state of expression (depending on gene function; dominant - recessive). DNA controls can be directly used.
4.3	Process controls	e.g. buffer used for extraction; a marker targeting the cytochrome oxidase gene as an internal amplification marker	YES	 a) Negative process control(s) b) Positive DNA control(s) that can be the control varieties c) Internal amplification control in case of a presence/absence marker

Chapter	Elements in a Standard characteristic- specific molecular marker protocol	Example	Essential information for harmoni- zation	Remark
5	Preparations	<i>e.g.</i> Sampling of seedlings 4 days old followed by DNA extraction using CTAB method	NO	Depending on the method used. Not in the Test Guideline. Detailed protocol(s) can be provided as an example in annex or available on request from the institute that developed the marker
6	Performance or Technique of the method	e.g. conventional PCR, TETRA-ARMS, qPCR, KASP, amplicon sequencing See TG/44/11/rev3 – Ad 51: ii DNA marker test add 6	YES	
6.1	Particular conditions	<i>e.g.</i> PCR protocol describing primer, enzyme, dNTP concentrations, PCR cycle scheme	NO	Depending on the method used. Not in the Test Guideline. Detailed protocol(s) can be provided as an example in annex or available on request from the institute that developed the marker
6.2	Particular hardware or infrastructure	<i>e.g.</i> machines, commercial kits, manufactures of components, lot numbers of chemicals	NO	Depending on the method used. Not in the Test Guideline. Detailed protocol(s) can be provided as an example in annex or available on request from the institute that developed the marker
7	Observations	<i>e.g.</i> Bands on agarose gel (conventional PCR), Ct values (qPCR) Variant call based on sequencing reads	NO	Depending on the method used. Not in the Test Guideline. Detailed protocol(s) can be provided as an example in annex or available on request from the institute that developed the marker
7.1	Validity of the results	e.g. for qPCR, Check for typical exponential amplification curves. Check if the controls are as expected (negative controls = no signal; positive controls = shows expected signals for all fluorophores).	YES	Depending on the method used.
8	Interpretation of the test results	See TG/44/11/rev3 – Ad 51: ii DNA marker test add 8	YES	Relation between alleles and expressions (with its notes)
9	Validation of the method,	This protocol was validated by a ring-test with different laboratories (<i>e.g.</i> Interlaboratory Comparative Test Report, INVITE 2024).	YES	
9.1	Contact Examination Office	<i>e.g.</i> Naktuinbouw	YES	Contact of the institute that developed this protocol, Name of the service.

VI. FOLLOW-UP SURVEY AFTER APPROVAL

34. Validation of the marker is not fixed as new genetics can arise from the market. This is a continuous evaluation process. Specificity should be re-assessed after validation acceptance using double testing at least during the first year with observation method.

35. After the first year of acceptance of the protocol, morphological checks on about 10% of the new varieties must be performed.

[Annex follows]

TWM/2/17

ANNEX

Guidelines for the validation of a new characteristic-specific molecular marker protocol as an alternative method for observation,

A collaborative work





Amandine LE VAN April 2024



TWM/2/17 Annex, page 2

Context

H2020 INVITE Project Task 5.2: Interlaboratory Validation of Tomato Resistance Markers

- Goal of this task: to validate the **reproducibility** of molecular markers linked to *Tomato mosaic virus* (ToMV), *Tomato spotted wilt virus* (TSWV) and *Fusarium oxysporum* f. sp. *lycopersici* Fol: 1EU/2US resistance/susceptibility in tomato (*Solanum lycopersicum*) and tomato rootstock.
 - Reproducibility in this study was evaluated as the consistency of results for each of the markers in the 3 participating laboratories CREA-DC, GEVES and Naktuinbouw
 - Specificity of the marker (correlation with the bioassay) previously evaluated in one lab (Naktuinbouw)
- The results were fully consistent within each lab and were 100% concordant for each of the markers in the three participating laboratories.

Groupe d'Étude et de contrôle des Variétés Et des Semences

Follow-up of this task:

- "Gold standard Protocol"
- Guideline on the validation of a new protocol based on molecular-marker



[End of Annex and of document]